

How to activate ZMAN license code

1. Download recent ZMAN software version and LabView runtime engine on <http://www.zivelab.com> or <http://www.wonatech.com>

Support – Software download menu in zivelab.com

ZIVE LAB

[About](#) [Site Map](#)

- ▶ Home
- ▶ Products
- ▶ **Support**
- ▶ News & Event
- ▶ Contact

Software Download



Support

- ▶ ZIVE Catalog
- ▶ WonATech Catalog
- ▶ Manual Download
- ▶ **Software Download**
- ▶ ScreenCasts
- ▶ FAQ
- ▶ Application/Tech Note
- ▶ Technical Support

User registration

*Institute:

Title:

Department:

*Name:

*Email:

*Country:

Phone:

FAX:

Model:
Select a Model... ▼

Serial No:

fields with an * are required

Software Download

marked link requires ID and password to download them.

If you are user and have a send serial number for ZIVELAB products, Please send the following "software request form" then we will promptly send software download FTP server's ID and password via email.

[Click here "Software Request Form"](#)

You can download other software by clicking following link.

Control/Data Acquisition Software for ZIVELAB products

- ZIVE SM software([ver6726](#)) for ZIVE SP/MP/BP/PP series with manual (60MB) [UPDATE](#)
- ZIVE Z100 Navigator software v.2.4 for Zcon/Z#/Z100 control (requires NI runtime engine 2013 32bit)
- ZIVE BM software([ver4121](#)) for BZA(Battery Impedance Analyser) series [UPDATE](#)

USB Driver / Installer

- * This driver is for ZIVE Z#,Zcon series
- ZIVE Device Centre ver1.5 Beta6
Do not use Device Center for SP1,SP2,SP5,MP1,MP2,MP5,PP1.

Data Analysis/tool Software

- **ZIVE ZMAN impedance data analysis software ([ver2.5](#))**
- ZIVE IVMAN DC data analysis software ver1.5 [UPDATE](#) Jan 2021
including:
 - ZIVE IVMAN Extractor module
 - ZIVE DA differential analysis software module
 - ZIVE PA Photovoltaic cell analysis software module
 - ZIVE PF peak find module
 - ZIVE TA Tafel analysis module
- ZIVE SIM4U digital CV simulation software (Freeware)[Oct 16, 2017](#)
- ZIVE Pathfinder software for EIS data file finding
- ZIVE Data Manager software ver1.28 [July 20, 2021](#) [UPDATE](#)
 - * Above softwares needs 32bit Labview runtime engine 2013 even if your OS is 64bit. 64bit Labview runtime is not available for these softwares.

LabVIEW RTE 2013 (32 bit): <http://www.ni.com/download/labview-run-time-engine-2013/4059/en/>

(copy the link and paste it into the URL bar.)

- * ZMAN/IVMAN/SIM4U software needs Microsoft .NET framework4.
<https://www.microsoft.com/en-us/download/details.aspx?id=17718>

Support

- Application Note
- Papers
- Software Download
- Manual Download
- Q&A
- WonATech Catalogs
- ZIVELAB Catalogs
- ScreenCasts

Software Upgrade

Support | Software Download

Software Download

marked link requires ID and password to download them.
 If you are user and have a send serial number for ZIVELAB products,
 Please send the following "software request form" then we will promptly send software download
 FTP server's ID and password via email.

[Click here "Software Request Form"](#)
 You can download other software by clicking following link.

Control/Data Acquisition Software for ZIVELAB products

-  [ZIVE SM software\(ver6726\)](#) for ZIVE SP/MP/BP/PP series with manual (60MB) [UPDATE](#)
-  [ZIVE Z100 Navigator software v.2.4](#) for Zcon/Z#/Z100 control (requires NI runtime engine 2013 32bit)
-  [ZIVE BM software\(ver4121\)](#) for BZA(Battery Impedance Analyser) series [UPDATE](#)

USB Driver / Installer

- * This driver is for ZIVE Z#,Zcon series
-  [ZIVE Device Centre ver1.5 Beta6](#)
Do not use Device Center for SP1,SP2,SP5,MP1,MP2,MP5,PP1.

Data Analysis/tool Software

- [ZIVE ZMAN impedance data analysis software \(ver2.5\)](#)
- [ZIVE IVMAN DC data analysis software ver1.5](#) [UPDATE](#) Jan 2021
including:
 - ZIVE IVMAN Extractor module
 - ZIVE DA differential analysis software module
 - ZIVE PA Photovoltaic cell analysis software module
 - ZIVE PF peak find module
 - ZIVE TA Tafel analysis module
- [ZIVE SIM4U digital CV simulation software \(Freeware\)](#) Oct 16, 2017
-  [ZIVE Pathfinder software for EIS data file finding](#)
- [ZIVE Data Manager software ver1.28](#) July 20, 2021 [UPDATE](#)


* Above softwares needs 32bit Labview runtime engine 2013 even if your OS is 64bit.
 64bit Labview runtime is not available for these softwares.

[LabVIEW RTE 2013 \(32 bit\): http://www.ni.com/download/labview-run-time-engine-2013/4059/en/](http://www.ni.com/download/labview-run-time-engine-2013/4059/en/)
 (copy the link and paste it into the URL bar.)


* ZMAN/IVMAN/SIM4U software needs Microsoft .NET framework4.
<https://www.microsoft.com/en-us/download/details.aspx?id=17718>

2. Install LabView runtime engine and ZMAN software on the PC which you want to use ZMAN. (Note: License code is per PC. You must install **the PC which you WANT to use ZMAN.**)
3. Send purchase order for ZMAN license to sales@wonatech.com

4. Run ZMAN software

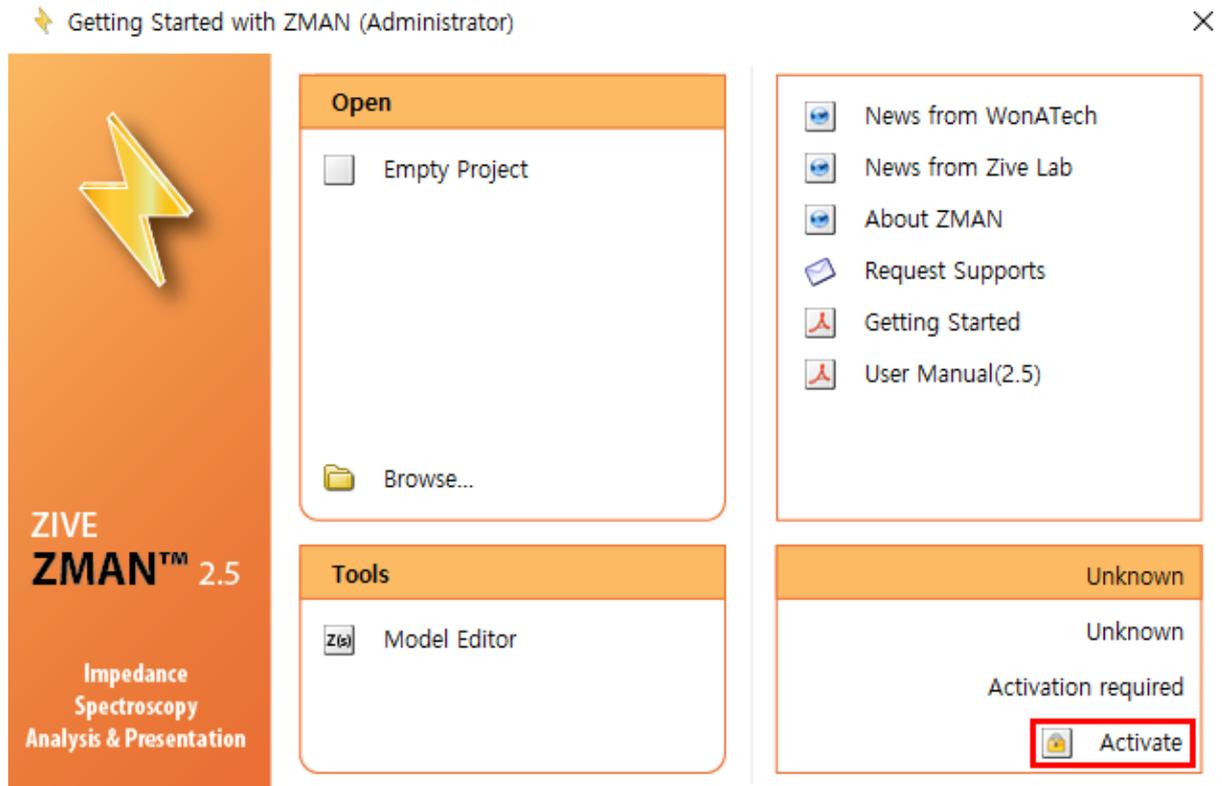

5. Click Activate

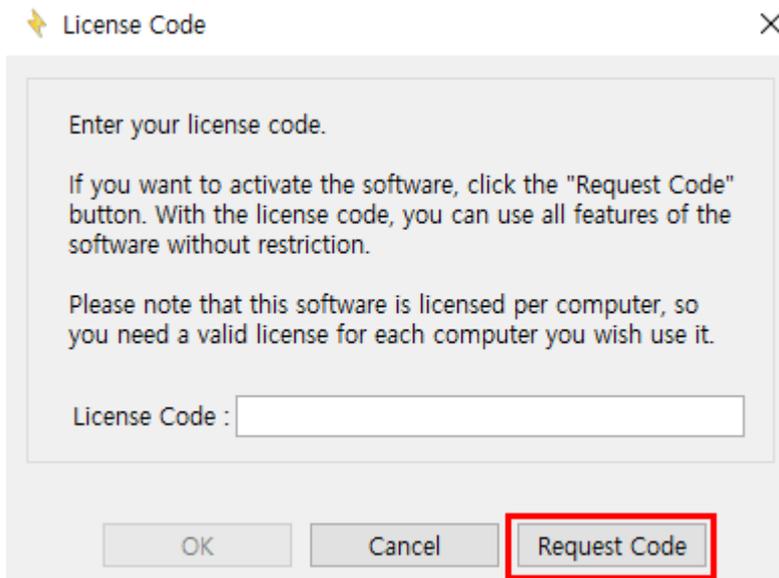

6. Click Request Code button


✦ User Info ✕

* Required


*First Name	*Last Name
<input type="text"/>	<input type="text"/>
*E-Mail Address	
<input type="text"/>	
*Phone Number	Fax Number
<input type="text"/>	<input type="text"/>
*Company/Institute	Department
<input type="text"/>	<input type="text"/>
*Street Address 1	
<input type="text"/>	
Street Address 2	
<input type="text"/>	
*Country	*Zip/Postal Code
<input type="text"/>	<input type="text"/>
Note	
<input type="text"/>	

Your right to privacy is a priority to us. We recognize your need for reasonable control of information that you share with us. Therefore, this statement is a corporate-wide guideline, and is not limited to our online business. We do not sell your information to outside organizations. Our goal in gathering information is to consistently improve our relationship with you, our customer.

7. Fill the information and click "OK" button

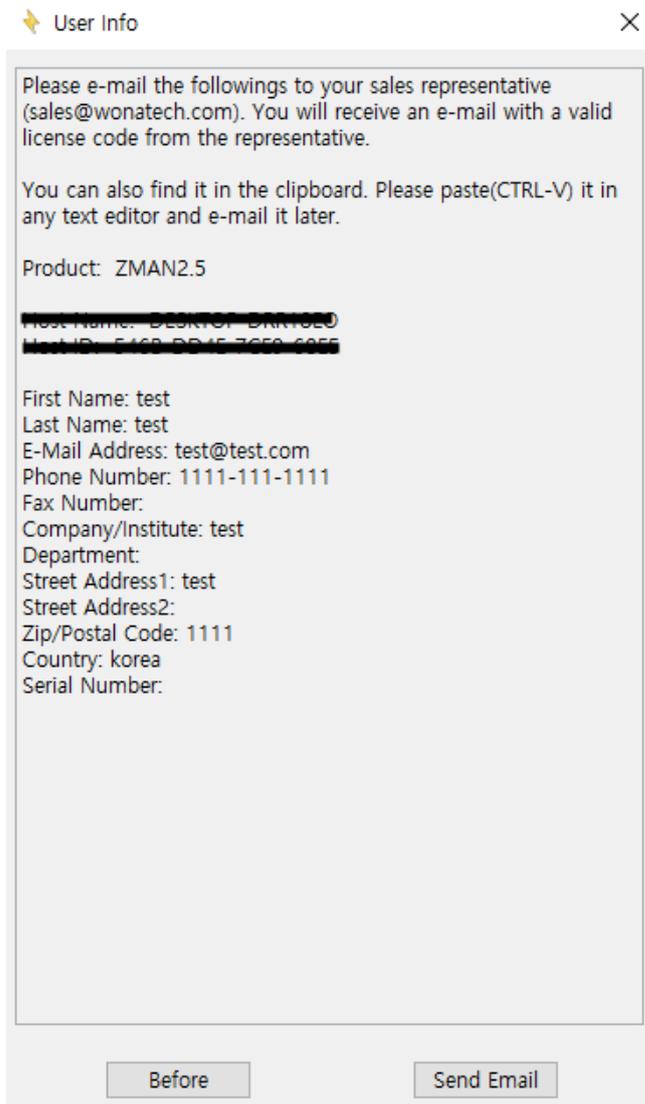


8. Click "send email" button or paste the contents using Ctrl+V on your email to sales@wonatech.com

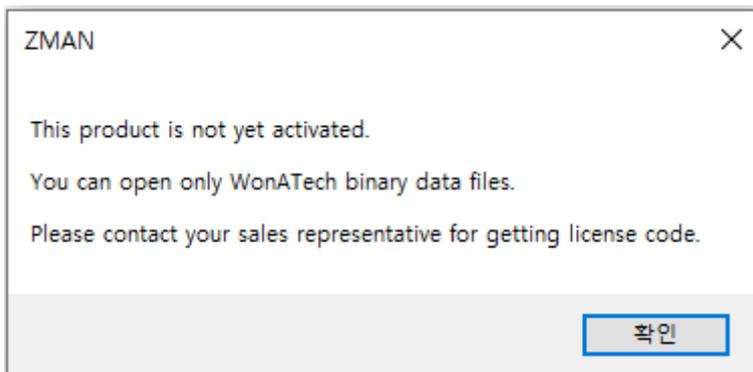


9. Wire transfer the order amount to WonATech account then we will send license code via email.
10. After receipt of license code, go to previous step 5 "Activate license code"


11. Input license code

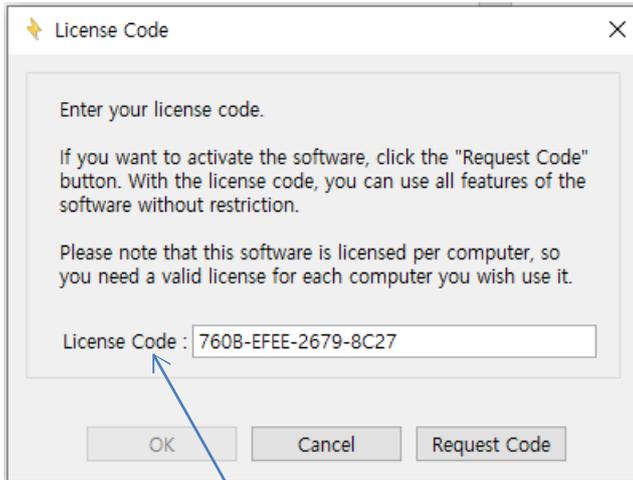

12. Click License code

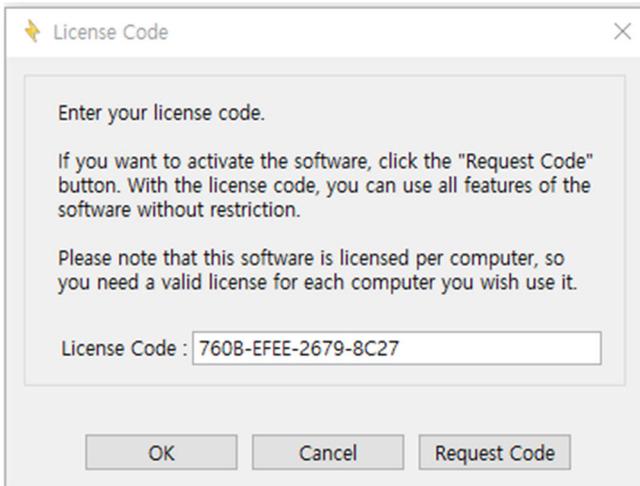

13. Click "OK button"

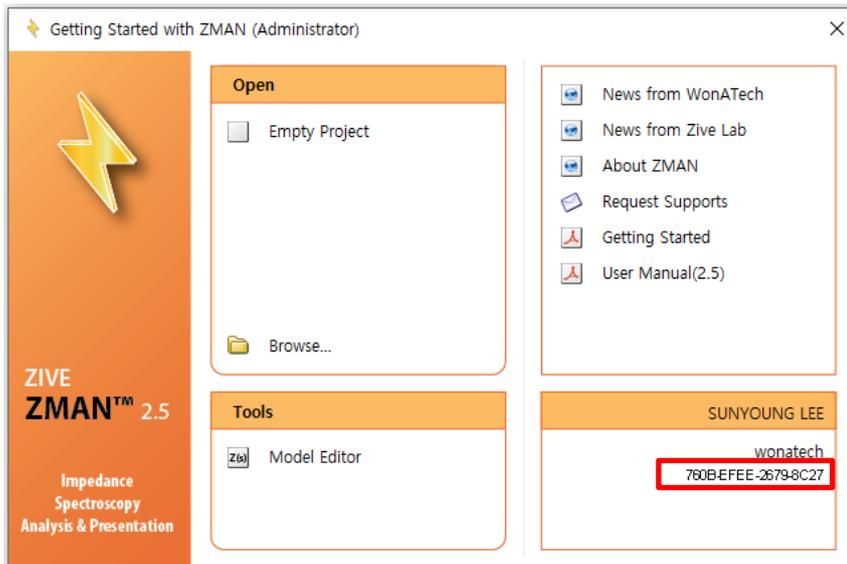

14. ZMAN is now activated.